

BIOCLEAN™

Laboratory Cleaning Solution



ISO 9001:2008
ISO 13485:2003

CE

BIOCLEAN™ is a detergent for Analytical, Pathological, R & D Laboratories where ultimate cleaning is essential for glass, plastic or other utensil.

USAGE

1. For floors, walls, benches, racks, shelves and lavatories, use a dilution of 100 mL per one liter of cold water and mop or brush for cleaning.
2. For cleaning glassware, bottles beakers, Porcelain dishes, use a dilution of 50 mL per one liter of warm water to get better result.
3. For microscopic slides, coverslip, wire meshes use a dilution of 100 mL per liter of hot water.

REAGENTS COMPOSITION

Akyle Sodium
benzene sulphinate 12%
Buffered to pH 7 ± 1

Revised 2.4.2014 KI

BIOCLEANENSURES

Best cleaning on any surface.

It is a stable, efficient surface active agent with a narrow range of pH variation. A 5% solution has a pH 6.5 to 7.5 that is near neutral pH.

Non-Corrosive, Non-toxic

PRECAUTIONS

BIOCLEAN is for laboratory use only.
Not for food, drug or household use.
Avoid contact with skin, eyes.
Keep away from children.

PACKING

1000 mL Plastic bottles and
5000 mL Plastic Jerry cans.

BIOLAB DIAGNOSTICS (I) PVT. LTD.

J-245, MIDC, Tarapur, Boisar – 401 501, MS.

E-mail : biolab@vsnl.com

www.biolabdiagnostics.com

Customer Care : (+ 9122) 28088243

BIOCLEAN™

Laboratory Cleaning Solution



BIOCLEAN™ is a detergent for Analytical, Pathological, R & D Laboratories where ultimate cleaning is essential for glass, plastic or other utensil.

BIOCLEAN ENSURES

Best cleaning on any surface. It is a stable, efficient surface active agent with a narrow range of pH variation. A 5% solution has a pH 6.5 to 7.5 that is near neutral pH.

one liter of cold water and mop or brush for cleaning.

2. For cleaning glassware, bottles beakers, Porcelain dishes, use a dilution of 50 mL per one liter of warm water to get better result.
3. For microscopic slides, coverslip, wire meshes use a dilution of 100 mL per liter of hot water.

REAGENTS COMPOSITION

Axyle Sodium

benzene sulphonate 12%

Buffered to pH 7 ± 1

BIOCLEAN is for laboratory use only.

Not for food, drug or household use.

Avoid contact with skin, eyes.

Keep away from children.

PACKING

1000 mL Plastic bottles and

5000 mL Plastic Jerry cans.

BIOLAB DIAGNOSTICS (I) PVT. LTD.

J-245, MIDC, Tarapur, Boisar – 401 501, MS.
E-mail : biolab@vsnl.com / www.biolabdiagnostics.com
Customer Care : (+ 9122) 28088243